## NOTICE D'INSTRUCTION - GANT DE BOITE A GANT EN BUTYL HAUTES PROPRIETES/EPDM TRITIUM (BHP/Eb TRI) 7/10mm MONTE SUR BAGUE SUPPORT

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## INSTRUCTIONS - GLOVEBOX GLOVES MADE OF BUTYL HIGH PROPERTIES/EPDM TRITIUM (BHP/Eb TRI) 7/10mm MOUNTED ON SUPPORT RING

### ➤ Scope/Characteristics of the material and expiry

- This glove can be used in nuclear industries, pharmaceutical industries and miscellaneous industries.
- This glove is manufactured using high-performance isobutylene polyisoprene (BHP) 80%, covered by EPDM (Ethylene-Propylene-Diene-Monomer (Eb)) 20%
- This glove protects against certain chemical and mechanical risks, against radioactive contamination and provides protection against micro-organisms (not viruses). It can be used in an atmosphere containing ozone. This glove does not protect against ionising radiation.
- The shelf life of this glove in its original packaging stored under the conditions described below is 3 years.

### ➤ Legislation, Risk analysis and recommendations for use

- This glove conforms to regulation 2016/425
- Please note, according to directive (UE) 2019/1832, the necessity to carry out a risk analysis relating to wearing the glove (e.g. breaking of leak tightness)
- Before each use, it is the responsibility of the user to check the quality and the integrity of the glove. In case of any doubt, it is recommended to discard the gloves
- For the assembly and disassembly of the gloves on the gloveboxes, please follow the instructions drafted by the safety manager and the glove port manufacturer.
- In case of a contamination, follow the instructions drafted by the safety manager.
- The gloves should be used at ambient temperature. Please contact the manufacturer for use under other temperature conditions.
- The gloves must be worn on clean and dry hands, with the nails cut short. Avoid wearing jewellery.
- It is recommended to not wear gloves when there is a risk of being caught by moving machine parts.
- During the use of unspecified chemical products (those not mentioned in the list of chemical products), please contact the manufacturer for more information.
- This glove can only be mounted if the client has a J2L system that allows receiving the glove (System not validated by IFTH during the EC type examination)

### ➤ Components:

- This glove does not contain substances in proportions in which they are known or suspected to have harmful effects on the user's health or hygiene under the foreseeable conditions of use.

### ➤ Properties of the glove

- *Chemical permeation EN ISO 374-1+A1 :2018*
- *resistance to deterioration EN374-4 :2013*

Tested products		Level of permeation (palm and cuff)	Average deterioration – palm (%)	Average deterioration – cuff (%)	Pictogram
K	Caustic soda 40%	6 of 6	-0.1	-8.8	
M	Nitric acid 65%	6 of 6	11.1	5.1	
P	Hydrogen peroxide 30%	6 of 6	6.6	8.6	

*This information does not reflect the actual duration of protection at the workplace, or the differentiation between the mixtures and pure chemical products. The chemical resistance has been evaluated under laboratory conditions using the samples collected in the palm and the cuff and only concerns the tested chemical product. It can be different if it is used in a mixture. It is recommended to check that the gloves are suitable for the intended use because the conditions at the place of work can differ from the typical test conditions, depending on the temperature, abrasion and deterioration. When they are worn, the protective gloves provide less resistance to hazardous chemical products owing to the change in their physical properties. Movements, tears, friction or deterioration caused by contact with chemical products, etc. can considerably reduce the actual duration of use. For corrosive chemical products, deterioration can be the most important factor to be taken into account in the selection of chemical-resistant gloves. Before use, it is recommended to inspect the gloves in order to ensure that they have no defects or imperfections.*

- *Protection against micro-organisms EN ISO 374-5: 2016*

Tests	Level	Pictogram
Airtightness	Compliant	
Water-tightness	Compliant	

*The resistance to penetration has been evaluated under laboratory conditions and only pertain to the tested test specimen.*

# INSTRUCTIONS - GLOVEBOX GLOVES MADE OF BUTYL HIGH PROPERTIES/EPDM BLACK TRITIUM (BHP/Eb TRI) 7/10mm MOUNTED ON SUPPORT RING

- Mechanical Strength **EN388+A1-2018**

Tests	Level	Pictogram
Abrasion	2 of 4	 2XX1X
Cut by slicing	X of 5	
Tear	X of 4	
Perforation	1 of 4	
X: test not carried out		

The material comprises two materials (two-coat). Hence, the overall classification does not necessarily reflect the performance of the outermost layer.

- Precision, Integrity and ozone cracking

Tests	Level
Precision <b>EN420+ A1:2009</b>	3 of 5
Integrity (pressure 30 mbar) <b>EN421-2010</b>	Compliant
Ozone cracking resistance <b>EN421-2010</b>	3 of 4

The highest class corresponds to the highest performance level

➤ **Integrity EN421-2010**

GLOVE PORT DIAMETER (mm)	FORM	ORIGINAL LEAK-TIGHTNESS PRESSURE (mbar)
136	54	20
156	50	30

The "glove port diameter" and "form" lists are not exhaustive. For any requests, please contact PIERCAN.

➤ **Marking**

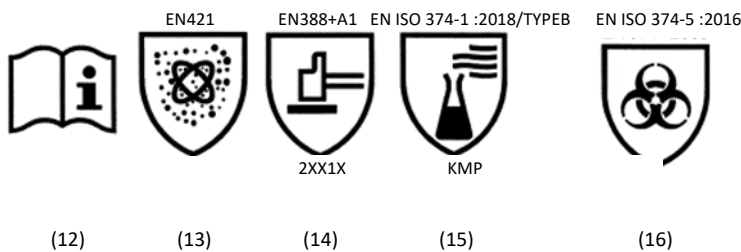
- Example of marking put on the glove and the packaging (see below)

PIERCAN LOGO

PIERCAN (1)

**0333** (2) **50750 BHP/Eb TRI 10 E7 GJABS [XXXXX]**  
 (3) (4) (5) (6) (7) (8) (9)  
**FR XXXXXX BATCH:XX XX** (10)  
**EXPIRY : XX/XX/XX** (11)

No.	DESCRIPTION
1	Manufacturer – In charge of release to the market
2	"CE" marking and No. of the body in charge of the annual inspection
3	Form
4	Length in mm
5	Material
6	Size
7	Thickness in 10th of mm
8	Glove mounted on ring
9	Product Item Code
10	Case number and batch number ("FR" French manufacturing "US" American manufacturing)
11	Date of expiry in storage
12	"INFORMATION" pictogram
13	"Protection against radioactive contamination" pictogram EN 421-2010
14	"Protection against mechanical risks" pictogram EN 388+A1 -2018
15	"Protection against chemical risks" pictogram EN ISO 374-1 +A1:2018
16	"Protection against micro-organisms" pictogram EN ISO 374-5 :2016

## INSTRUCTIONS - GLOVEBOX GLOVES MADE OF BUTYL HIGH PROPERTIES/EPDM BLACK TRITIUM (BHP/Eb TRI) 7/10mm MOUNTED ON SUPPORT RING

➤ **Size available in the main ambidextrous model**

FORM	∅ RDG (glove port diameter)	∅ BS (internal diameter of the support ring)	SIZE					
			7.5	8.5	9	9.5	10	10.5
50 (GJABS)	136	131					x	

➤ **Suitable packaging for transport**

- The gloves should be transported in their original packaging.

➤ **Storage**

- Store in a dry place, away from light
- Storage temperature: 5 to 25°C.
- Do not store near electrical installations to avoid accelerated ageing

➤ **Cleaning/decontamination and maintenance**

- If necessary, the gloves are washed with soapy water, rinsed with clear water, and dried under a low temperature air flow (<40°C) and powdered. The performance is not guaranteed by the IFTH after washing
- Do not use chemical products
- If the use of unspecified chemical products is necessary (those not mentioned in the list of chemical products), please contact the manufacturer for more information.
- Do not use pointed or sharp objects like wire brushes, sandpaper or similar objects

➤ **Treatment of the glove**

- Non-recyclable gloves
- It can be discarded as simple non-toxic waste if it is not soiled by a hazardous product
- If a glove is soiled, discard it through the appropriate channel

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